



STERILIZER (STEAM ONLY)	MAX TEMPERATURE	MAX PRESSURE	MAX TIME
Steam autoclave	136 C (276 F)	15psi	15min

- Laboratory Grade Hand Crimpers and Decrimpers – Remove Grips prior to process
- Painted Hand Crimpers and Decrimpers- Not Recommended
- Power Crimper, Heads, and stand only (Do not autoclave regulators, hoses, foot pedals, verification gauge)
- VAC Units and Heads only (Do not autoclave regulators, hoses)

There will be a slight discoloring of the surface areas and the lifetime of all units will be decreased. Autoclaving will void warranties.

Kebby Products have not been evaluated for clean room classification, and therefore, no cleanroom classification has been assigned. It is up to the management of the cleanroom facilities or clean areas to determine if Kebby Products are suitable for their applications.